

## SONICA<sup>®</sup> 2400 S3

Product name	Code
<b>SONICA 2400 M S3 *</b>	090.003.0006
<b>SONICA 2400 MH S3 *</b>	090.003.0007
<b>SONICA 2400 ETH S3 *</b>	090.003.0008
<b>SONICA 2400 EP S3 *</b>	090.003.0009
<b>Ultrasonic Cleaner 4,5 Litres</b>	

### Accessories

Stainless steel wire basket Internal dim. 265x120x45mm	090.004.0003
Stainless steel lid 2400	090.004.0019
Stainless steel auxiliary tank SONICA 2400 Indicative dim. 300x150x100mm	090.004.0041
Stainless steel beaker holder for two glass or plastic beaker	090.004.0036
Plastic beaker 250 ml	090.004.0032
Glass beaker 400 ml	090.004.0033

### Further accessories on request



\* *Stainless steel tank made in AISI 304 thickness 08/10 with round tank corners for a better removing of dirt residuals after cleaning process. The tank is conical and realized with a press machine*

### General Characteristics

Ultrasonic Cleaning is a safe and efficient modern procedure, which ensures perfect cleaning within the shortest time possible. It has many advantages over ordinary cleaning methods. Ultrasonics can remove even the most tenacious deposits from parts which would be difficult to reach by hand, it also reduces the risk of injury or infection which may occur during manual cleaning.

### Notes for use as "Medical Devices"

**Denomination:** Ultrasonic Cleaning Device UMDNS-nomenclature (ECRI) 14-263

**Field of application:** Cleaning and disinfection of surgical and dental instruments

**Classification** (Medical Devices Directive 93/42/EEC, appendix IX): Class I, Non Active MD, Non Invasive MD, Non Implantable MD

### "Medical Device" registration and classification

**Registration Number:** 103291/R SOLTEC Srl **CND Classification:** Z12011302 Ultrasonic Cleaning Device

#### SOLTEC S.r.l.

Uffici Commerciali e Sede Legale: Via G. Röntgen 16 - 20136 Milano  
 Tel. +39 02 58 308 378 - Fax +39 02 58 308 595 - [www.soltec.it](http://www.soltec.it) - [info@soltec.it](mailto:info@soltec.it)  
 Produzione e Magazzino: Via G. Castelbarco, 17 - 20136 Milano - Tel. +39 02 58 324 131

Cod. fisc. e P.IVA: IT11127210158  
 N. Mecc.: MI 221750 - C.C.I.A.A. Milano N. 1441548  
 Iscr. Trib. Mi Reg. Soc. N. 343391/8444/41  
 Cap. Soc. Euro 110.500 i.v. - Iscr. Registro AEE Nr. IT08020000000435



Per la gestione dei RAEE  
 questa azienda si affida a  
**Consorzio ecoR'it**



ISO 9001:2015  
 Reg. N.: 9319-A  
 ISO 13485:2016  
 Reg. N.: 9319-M

# SONICA<sup>®</sup> 2400 S3

Technical specification subject to change

## Technical Data

Main line voltage	230/240V~50/60Hz
Main input power/consumption	130 W
Main input power with heating	305 W
Peak HF Ultrasound Power	260 W
US frequency	39 kHz ± 1 kHz with SWEEP SYSTEM Technology
Liquid drain	tube 10x8mm
Transducers	2 (2 Piezoelectric ceramics each)
Tank capacity [L]	4,5
Weight [KG]	3,8
Installation category	Class II according to EN 61010-1
Environmental conditions	Temperature Rel. Hum. 80% up to 31°C with linear decrease up to 50% at 40°C
IP grade	IP 22
CE-compliant	2004/108/EC EMC Directive (EN61326) 2006/95/EC LV Directive (EN 61010-1) 93/42/CEE Medical Device Directive Risk Class I according to the rule 12 to the Directive 93/42/EEC Annex IX

## Approximately Internal Tank Dimensions


W/D/H (mm)	300 / 150 / 100
Max filling volume tank	4,5 L

## External Appliance Dimensions

W/D/H (mm)	325 / 175 / 260
Max operating tank volume	3,15 L

## Constructive Materials

Housing	Stainless steel AISI 304
Tank	Stainless steel AISI 304

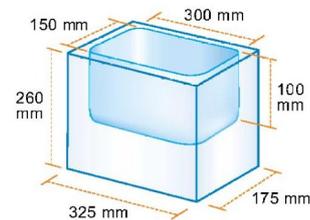

## Control panel available series



**M serie:** The simplest version in the SONICA range is equipped with a timer which can be set for 15 minutes. It can also be set to manual operation up to 1 hour of continuous. No heating system  
**Sweep System** generator

**MH serie:** This version is equipped with a timer which can be set for 15 minutes. It can also be set to manual operation up to 1 hour of continuous. The thermostat for the heater is set at 60°C.  
**Sweep System** generator

**ETH serie:** Cleaning times and temperature controlled electronically by a microprocessor. Ultrasonic cleaning times 5-10-15 minutes. Double washing times 10-20-30 minutes by pressing twice the minutes key  
**Sweep System** generator

**EP serie:** The most technologically advanced of the SONICA range: **Sweep System** generator; modern digital display control panel window; degas time setting; cleaning times from 0 to 99 minutes; tank liquid temperature setting up to 70°C; memory for 9 cleaning programmes; jet program function to start quickly and easily cleaning cycle.